


# Dispenser arm kit for Frac-950

## Instructions

### Removing the standard mode components

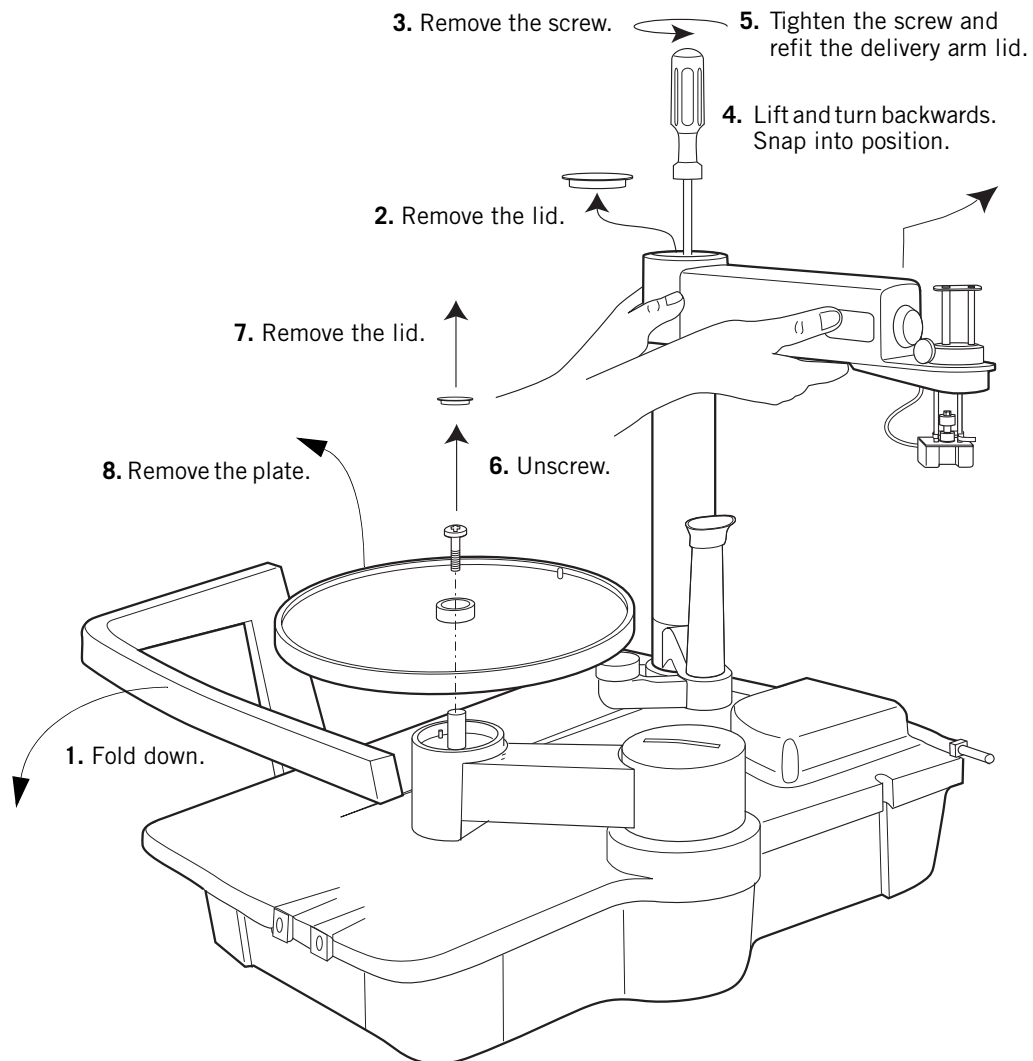


### Installing the rotating shaft

2. Use the keying guide to position the shaft correctly.

3. Insert and tighten the screw.

1. Fit the shaft to the guide pin.

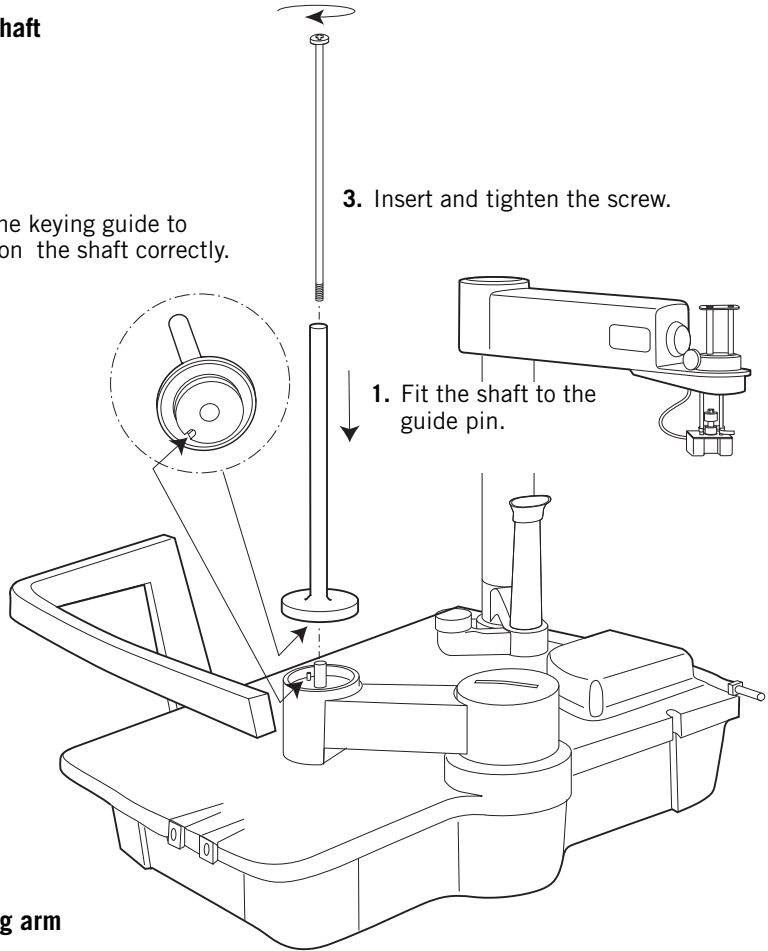


### Mounting the dispensing arm

1. Loosen the knurled screw.

2. Fit the tip of the knurled screw to the shaft recess.

3. Tighten the knurled

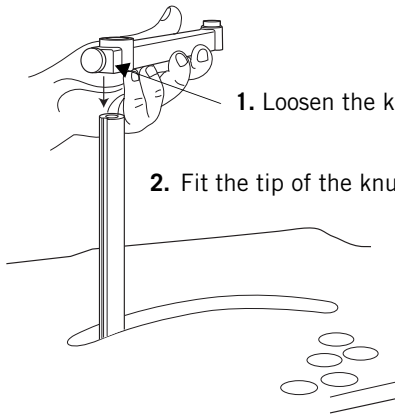


## Installing the capillary

3. Thread the free end of the capillary through the fingertight connector and the swivel. Allow the capillary end to stick out 1-2 mm. Hold the swivel, with the 8 mm key supplied, and tighten the connector.

2. Fit a fingertight connector to the swivel.

1. Connect the capillary to the accumulator.

4. Adjust the height of the vessels (a distance between the arm and the vessels of 0.5-1 cm is appropriate). Tighten the knurled screw.

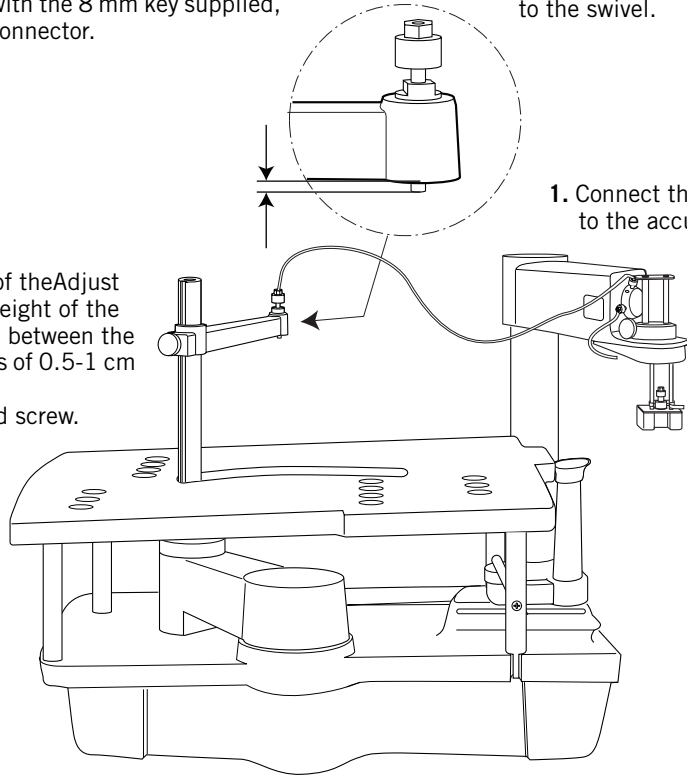

## Important user information

### Declaration of conformity

The **CE** symbol and corresponding declaration of conformity is valid for the instrument when it is:

- connected to other CE-marked Amersham Biosciences instruments, or
- connected to other products recommended or described in this manual, and
- used in the same state as it was delivered from Amersham Biosciences except for alterations described in this manual.

### Terms and Conditions of Sale

Goods and services are sold subject to the terms and conditions of sale of the company within the Amersham Biosciences group which supplies them. A copy of these terms and conditions is available on request.

Should you have any comments on this product, we will be pleased to receive them at:

**Amersham Biosciences AB**  
SE-751 84 Uppsala  
Sweden

### Trademarks

Amersham and Amersham Biosciences are trademarks of Amersham plc.

### Office Addresses

**Amersham Biosciences AB**  
SE-751 84 Uppsala  
Sweden

**Amersham Biosciences UK Limited**  
Amersham Place, Little Chalfont, Bucks, HP7 9NA  
England

**Amersham Biosciences Corp.**  
800 Centennial Avenue  
PO Box 1327  
Piscataway, NJ 08855  
USA

**Amersham Biosciences Europe GmbH**  
Munzinger Strasse 9  
D-79111 Freiburg  
Germany

**Amersham Biosciences KK**  
Sanken Building  
3-25-1 Hyakunincho, Shinjuku-ku  
Tokyo 169-0073  
Japan

© Amersham Biosciences AB 2002  
– All rights reserved