

Confocal Application Notes

Vol. 4 March 2006



CROSSTALK REMOVAL

Prepared by Myriam Gastard, Ph.D.
Exton Support Group, Leica Microsystems, Inc.



One of the most frequently asked questions from our confocal users is how to correct for crosstalk during simultaneous image acquisition. A precondition to correct for crosstalk is a confocal microscope equipped with an AOTF (Acousto Optical Tunable Filter) and a spectral detection system like in the SP5.

The AOTF is necessary to attenuate every laser line individually from 0%-100% and the spectral detection system enables you to create your own detection filter settings optimized for crosstalk removal.

This application note will demonstrate how to reduce or eliminate crosstalk using the two fluorophores FITC and MitoTracker Red. This protocol can be used to help eliminate crosstalk between virtually any two or three fluorophores. With the Leica SP5 system these settings can be saved for future use to optimally separate the two fluorophores.

PROTOCOL:

1. To verify whether you do have crosstalk or not in a double or triple labeled specimen e.g. using FITC/Mito Red you should begin with the simultaneous detection of both labeling. (Emission/Excitation of FITC: 494/518 nm and Mito Red: 578/599 nm).

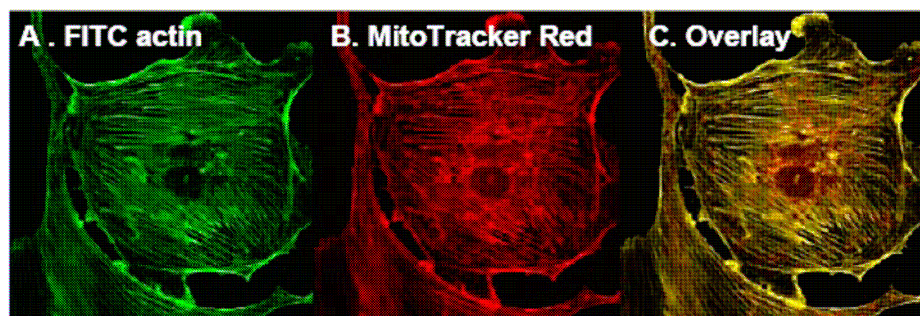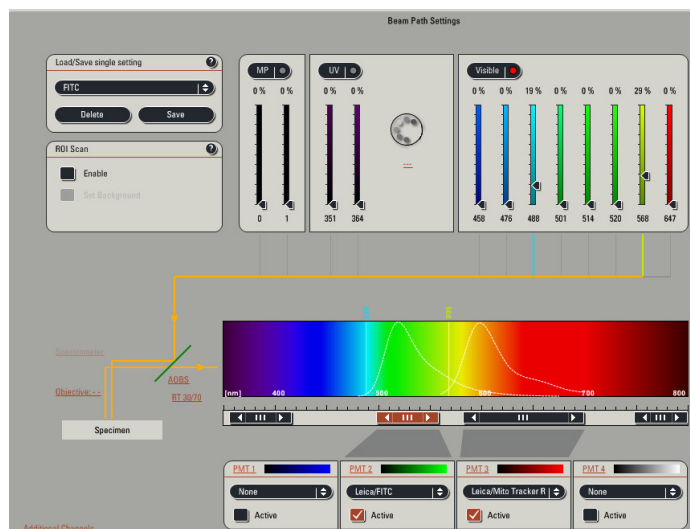

Leica Microsystems Inc.
410 Eagleview Blvd, Ste. 107
Exton, PA 19341



Telephone (610)-321-0460
Toll Free 866-830-0735
Fax (610)-321-0425
www.confocal-microscopy.com

Confocal Application Notes

Vol. 4 March 2006



2. Now reduce the excitation light source of the MitoTracker Red label (red channel) down to 0%. Any structure that you still see in the red channel does not arise from the MitoTracker Red label but is the red fraction of the fluorescein emission and can be categorized as crosstalk.

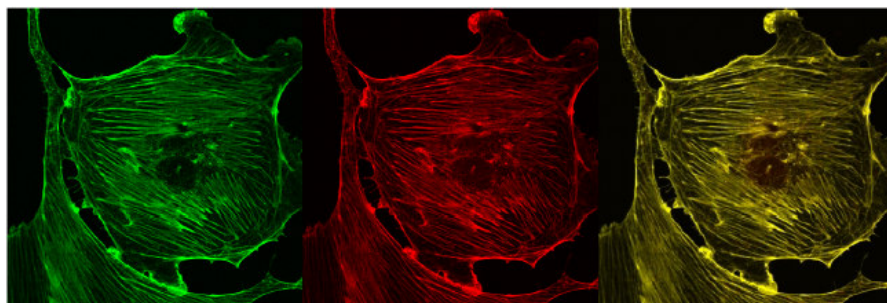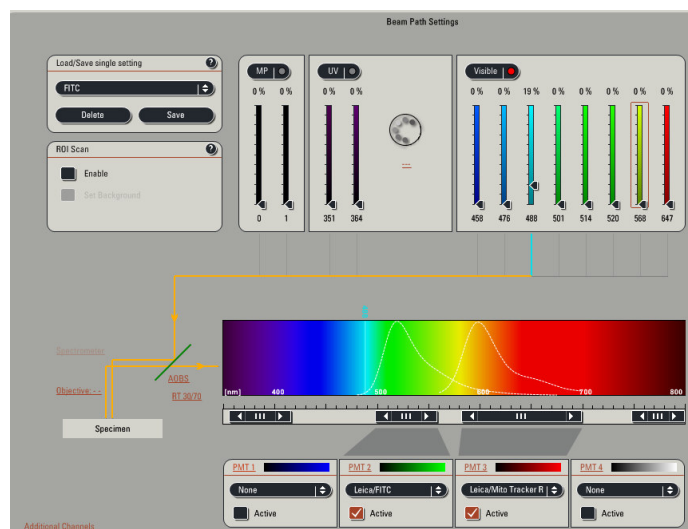

(Crosstalk from FITC into the MitoTracker Red channel)

Leica Microsystems Inc.
410 Eagleview Blvd, Ste. 107
Exton, PA 19341


Telephone (610)-321-0460
Toll Free 866-830-0735
Fax (610)-321-0425
www.confocal-microscopy.com

Confocal Application Notes

Vol. 4 March 2006



3. To eliminate these effects (i.e. the crosstalk) you must reduce the intensity of excitation light for FITC (488 nm) **and** move the detection slits of the MitoTracker Red detector (PMT3) away from the emission tail of FITC until the red signal disappears. Minor adjustments to the gain/offset of FITC may need to be made. If these steps don't resolve the crosstalk, then a sequential scanning can be setup to eliminate it (Will be addressed in our April Application Notes).

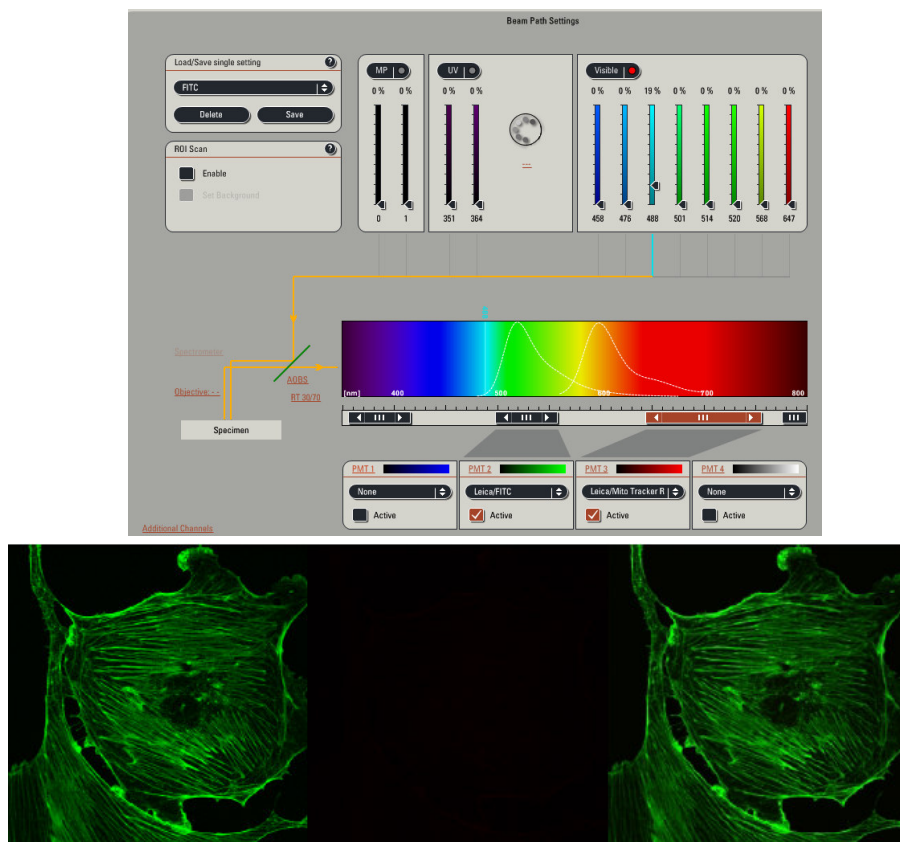

Leica Microsystems Inc.
410 Eagleview Blvd, Ste. 107
Exton, PA 19341


Telephone (610)-321-0460
Toll Free 866-830-0735
Fax (610)-321-0425
www.confocal-microscopy.com

Confocal Application Notes

Vol. 4 March 2006



4. Once you have found a satisfactory setup of the excitation light for FITC 488 nm laser line (i.e. good or acceptable signal-to-noise ratio for the FITC signal and little or no signal remaining in the red) you can now again excite MitoTracker Red with the 568 nm line. Adjust the intensity of the 568 nm line until you have a good signal-to-noise ratio.

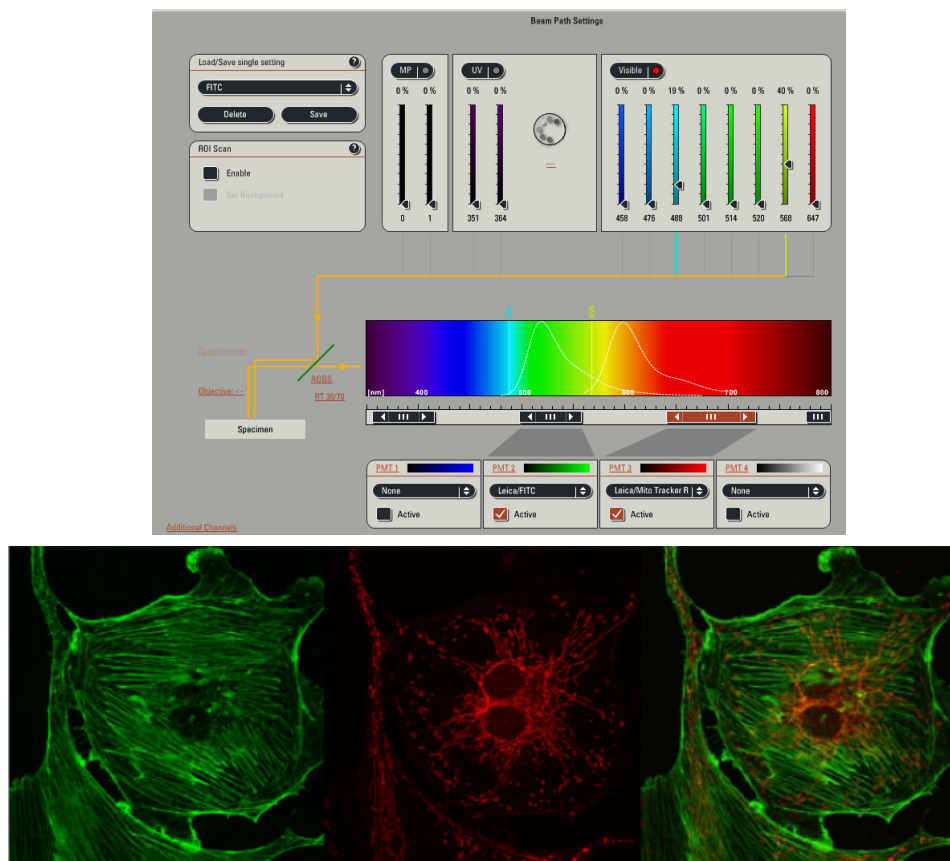

5. These settings can now be saved for future use when using these fluorophores.



Leica Microsystems Inc.
410 Eagleview Blvd, Ste. 107
Exton, PA 19341

Telephone (610)-321-0460
Toll Free 866-830-0735
Fax (610)-321-0425
www.confocal-microscopy.com